

1. Equipment Installation

STEP 1 Plug the fan heater cables into the respective connectors as per the depicted color code.

STEP 2 Connect the heating box cables to the Heating Box 1 and Heating Box 2 ports on the rear panel of the controller and enclosure.

#2 H203-HEATING-BOX-CABLE

STEP 4 Connect the OKO-TOUCH to the H203-T-UNIT by using the OKO-TOUCH-CABLE.

STEP 5 Pass the on-stage chamber connector through the cable-pass slot indicated and plug it to the Chamber port on the rear panel of the H203-T-UNIT.

STEP 13 Plug the controller into a power point.

Power Supply-Temperature Control Unit

STEP 15 Plug the enclosure into a power point.

ENCLOSURE POWER CORD

2. System Start & Quick Usage

TURN THE SYSTEM ON

Turn the system ON by pressing and holding the power button on the OKO-TOUCH.

SET DATE&TIME

Navigate to 'Settings' > 'System' > and click on the 'Date & Time' icon .

SETPOINT

Insert the values setpoint by clicking on the corresponding badge.

ALARMS

To set an alarm select the value of Deviation (tolerance) from the set point and the time for which the value must be out of tolerance to send out an alarm.

OVERVIEW PAGE

System Overview