

1. Equipment Installation

STEP 1 Connect the OKO-TOUCH to the H303-T-UNIT by using the OKO-TOUCH-CABLE.

OKO-TOUCH-CABLE

STEP 2 Install the chamber on the stage.

STEP 3 Plug the on-stage chamber connector to the Chamber port on the rear panel of the H303-T-UNIT-BL.

Chamber

STEP 4 Connect the TC-F to read the temperature inside the sample.

TC-F
TC-EXTENSION CABLE

Connect the TC to read the room temperature.

TC

STEP 5 Plug the controllers into a power point.

Power Supply-Temperature Control Unit

2. System Start & Quick Usage

TURN THE SYSTEM ON

Turn the system ON by pressing and holding the power button on the OKO-TOUCH.

SET DATE&TIME

Navigate to 'Settings' > 'System' > and click on the 'Date & Time' icon .

SETPOINT

Insert the values setpoint by clicking on the corresponding badge.

T set point 37.0°C

ALARMS

To set an alarm select the value of Deviation (tolerance) from the set point and the time for which the value must be out of tolerance to send out an alarm.

Settings Alarm Temperature

OVERVIEW PAGE

System Overview

Room 23°C

Lid 39.3°C

Base 37.1°C

ObH 36.6°C

T-Sensor 37.1°C